

PAWKIT



TABLE OF CONTENTS

1. Introduction.....	1
2. Operation	2
2.1 Setup	2
2.1.1 Sample Preparation	3
2.1.2 Taking a Measurement	4
3. System.....	9
3.1 Specifications.....	9
3.2 PAWKIT Components.....	10
3.3 Theory	11
3.3.1 Moisture Content	11
3.3.2 Water Activity	12
3.3.3 Water Potential	15
3.3.4 Sorption Isotherms	15
4. Service.....	16
4.1 Verification and Calibration	16
4.1.1 Verification Standards	16
4.1.2 Verify Calibration	17
4.2 Maintenance	19
4.2.1 Cleaning.....	19
4.2.2 Filter Replacement	21
4.2.3 Battery Replacement	22
4.3 Repairs	25
4.4 Troubleshooting	25

4.5 Customer Support.....	26
4.6 Terms and Conditions	27

References	28
-------------------------	-----------

Index	29
--------------------	-----------

1. INTRODUCTION

Thank you for choosing the AQUALAB PAWKIT Water Activity Meter.

The PAWKIT is designed to be a simple, rapid, and portable system for water activity measurement to ensure product safety. It is easy to use, durable, and requires little maintenance.

This manual includes instructions for setting up, verifying the calibration, preparing samples, and maintaining and caring for the PAWKIT. Please read these instructions before operating the PAWKIT to ensure that the instrument performs to its full potential.

Verify that all PAWKIT components are included and appear in good condition:

- PAWKIT main unit
- Durable carrying case
- Calibration certificate

2. OPERATION

Please read all instructions before operating the PAWKIT to ensure it performs to its full potential.

PRECAUTIONS

AQUALAB products are built to the highest standards, but misuse may damage the product and possibly void the manufacturer's warranty. Follow the recommended setup instructions and arrange proper protections to safeguard products from damage.

2.1 SETUP

Follow the steps listed in [Table 1](#) to set up the PAWKIT and start collecting data.

SAMPLING PRECAUTIONS

Sensor calibration can shift when the PAWKIT is exposed to a variety of volatile substances or to samples with water activities near 1.00 over a long period of time. To prevent a sensor calibration shift, always remove samples as soon as the PAWKIT is finished sampling (a beep signals when it is finished) to avoid damage to the sensor. If a sample is accidentally left in the chamber for an extended period of time, be sure to check the calibration when the instrument is used next.

Table 1 Preparation

Tools Needed	PAWKIT unit Material to sample Plastic sample cup WARNING: The PAWKIT is not intended for outdoor use.
Preparation	Select Clean, Level Location Always place the PAWKIT on a level surface when measuring to ensure correct and consistent operation. This also avoids the risk of sample material spilling inside the instrument and causing contamination. Place the PAWKIT in a location where the temperature remains fairly stable, to avoid inaccurate readings (away from air conditioner and heater vents, windows, etc.) IMPORTANT: Temperature fluctuations or vibrations will strongly affect the accuracy of measurement results. Sample Preparation Prepare the sample as described in Section 2.1.1 .

Table 1 Preparation (continued)

Installation	Sample Placement Place the PAWKIT over the sample cup as described
---------------------	--

2.1.1 SAMPLE PREPARATION

Carefully prepare samples to get the best readings possible. Always use clean sampling practices to ensure accurate and repeatable readings. Follow the steps below when preparing samples.

1. Make sure the sample to be measured is homogenous.

Multicomponent samples (e.g., muffins with raisins) or samples that have outside coatings (like deep-fried, breaded foods) can be measured, but may take longer to equilibrate. Samples like these may require additional preparation (cutting, crushing, or grinding) to obtain a representative sample.

2. Wash hands thoroughly or put gloves on prior to handling the PAWKIT and sample cup.
3. Fill the sample cup no more than half full.

Completely cover the bottom of the cup with the sample, if possible. The PAWKIT is able to accurately measure a sample that leaves small spaces of the cup bottom exposed. For example, raisins only need to be placed in the cup and not flattened to cover the bottom. A larger sample surface area increases instrument efficiency by shortening the time needed to reach vapor equilibrium. However, if the sample cup is too full, contamination of the sensor becomes a risk, which can lead to inaccurate readings.

4. Clean the bottom, edges, and rim of the sample cup thoroughly with KIMWIPES® strips ([Section 4.2](#)).

Wipe any excess sample material from the rim of the cup with a clean KIMWIPES tissue. Material left on the rim or the outside of the cup can be transferred to subsequent samples and may affect the reading accuracy. The rim of the cup forms a vapor seal with the sensor. Any sample material left on the cup rim may prevent this seal and contaminate future samples.

If a sample reading will be taken at a later time, put the sample cup disposable lid on the cup to restrict water transfer. To seal the lid, wrap tape or Parafilm® wrapper completely around the cup/lid junction. It is necessary to seal the cup if it is going to be a long time before measurements will be made again.

The PAWKIT makes its most accurate measurements when the temperature of the sample and instrument are within 1 °C. If the sample is too warm, the thermometer icon on the left of the screen appears ([Figure 1](#)).

OPERATION

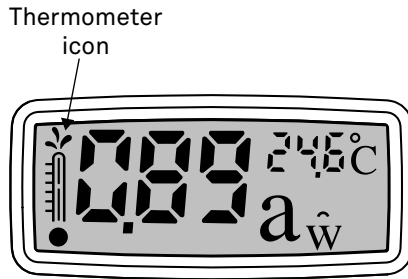


Figure 1 Thermometer icon on screen

The instrument beeps, when the sample temperature is too high, indicating there is a danger of condensing water in the sample chamber and on the sensor. If this warning appears while sampling, remove the PAWKIT, place the cup lid on the sample, and wait until the sample has reached ambient temperature before attempting to read again.

If the sample is colder than the ambient temperature of the PAWKIT, the accuracy of the reading after 5 min may be questionable. Wait until the sample temperature is similar to that of the PAWKIT.

2.1.2 TAKING A MEASUREMENT

1. Open the PAWKIT by holding the case near the LCD with one hand and pulling down on the plastic sensor cover tab with the other hand.

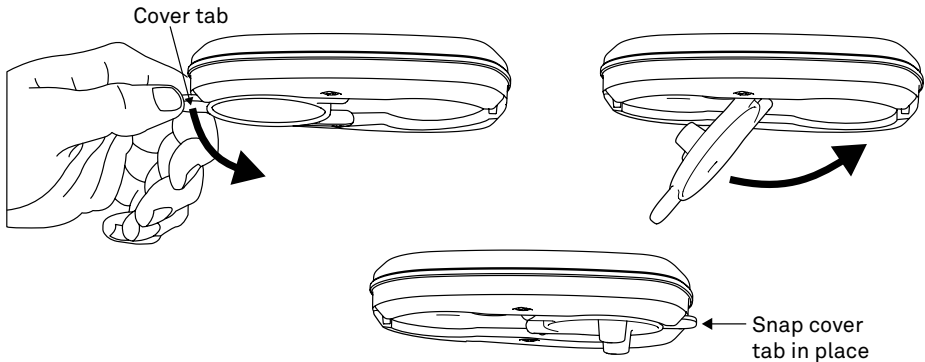


Figure 2 Open the PAWKIT

The sensor cover rotates and snaps into the open position.

OPERATION

2. Place the prepared sample cup onto a level surface.

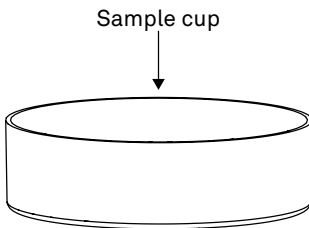


Figure 3 Sample cup placement

3. Place the opened PAWKIT onto the prepared sample cup.
The cup fits under the sensor into a recess in the bottom of the PAWKIT.

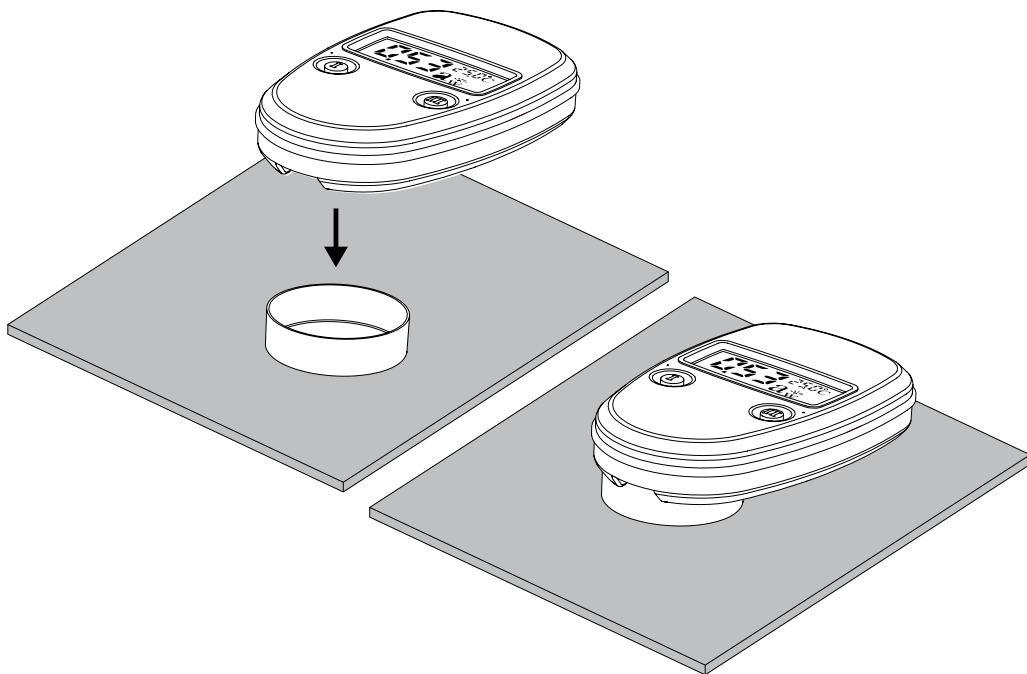


Figure 4 PAWKIT placement on sample cup

A correctly positioned cup results in the PAWKIT being level on the bench when sitting on the cup and the sensor cover legs. Ensure the rim of the cup is entirely within the recess of the sample chamber (Figure 4). Otherwise, the PAWKIT may not be level on the bench, and the cup might not make a vapor seal with the sensor.

OPERATION

4. Press the I button (located on the left) to turn on the instrument.

The instrument will display the last reading taken. This allows the PAWKIT to take a measurement without anyone attending the instrument throughout the measurement. If the PAWKIT is already on, proceed to the next step.



Figure 5 I button


5. Press the I button a second time to begin the water activity measurement. The LCD display resets to 0.00 a_w .

NOTE: Pressing the I button any time during a measurement restarts the water activity measurement.

Once the measurement process has been started, the PAWKIT begins to display water activity measurements as well as temperature after 5 s, updating the display every second thereafter.

During this time the sunburst icon to the right of the water activity value indicates it is measuring. As it measures, the beams of the sunburst move from left to right.

NOTE: The final water activity measurement is not displayed until the instrument beeps and the sunburst icon disappears from the screen.

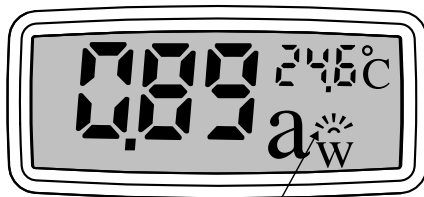

Sunburst and
beams icon

Figure 6 Sunburst and beams icon

OPERATION

If an error code of 9.99 is received at any time during the process, it indicates that the sensor has failed and that the instrument needs to be serviced. Refer to [Section 4.4](#) for instructions on how to return the PAWKIT for repair.

NOTE: DO NOT lift or move the instrument during the measurement because it could contaminate the chamber, break the chamber vapor seal, and invalidate the water activity measurement.

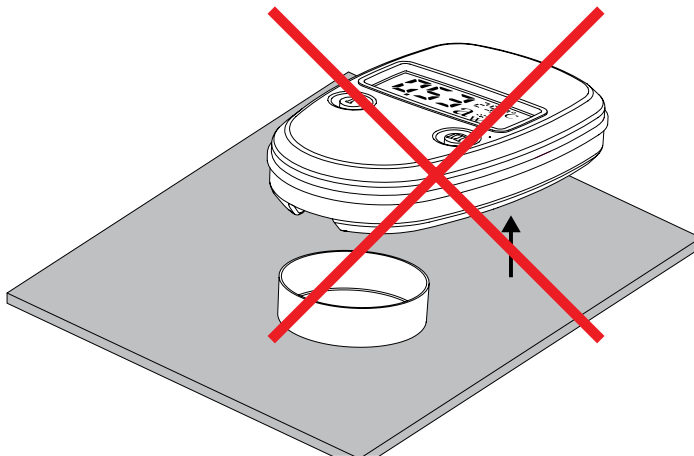


Figure 7 DO NOT lift or move instrument during measurement

6. After 5 min, the instrument displays the final water activity and beeps 5 times. The sunburst disappears when the water activity reading is finished.

NOTE: AQUALAB recommends recording the reading value before proceeding.

At this point, either restart the measurement by pressing **I** button again or end the measurement procedure.

7. Remove the sample cup by lifting the PAWKIT

Lift the PAWKIT straight up as shown in [Figure 8](#) to avoid spilling the sample cup. The sample may now be discarded or covered with a lid if it is to be remeasured at a later time.

OPERATION

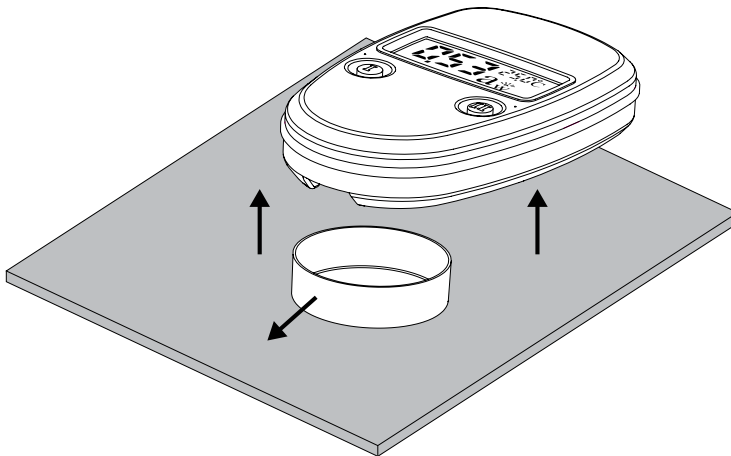


Figure 8 Remove PAWKIT when measurement complete

8. When finished taking readings, close the instrument.

With one hand holding the case near the LCD, pull down on the plastic sensor cover tab with the other hand and rotate until it snaps into the closed position covering the sensors.

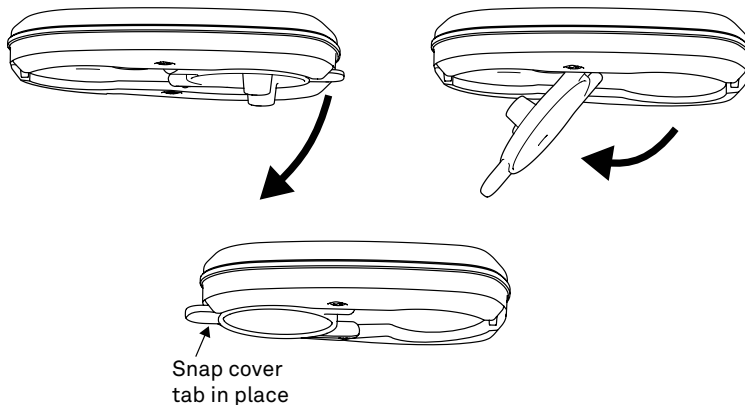


Figure 9 Closing the PAWKIT

To turn off the PAWKIT, leave it idle for more than 5 min and it will shut off automatically. If the PAWKIT has automatically shut itself off, pressing the I button wakes up the instrument and displays the last water activity measurement.

NOTE: Remember to close the lid before storing the case.

3. SYSTEM

This section describes the specifications, components, and theory of the PAWKIT.

3.1 SPECIFICATIONS

MEASUREMENT SPECIFICATIONS

Water Activity

Range	0.00-1.00 a_w
-------	-----------------

Resolution	0.01
------------	------

Accuracy	± 0.02
----------	------------

Sample Temperature

Resolution	0.1 °C
------------	--------

Accuracy	± 0.2 °C
----------	--------------

Read Time

5 min

PHYSICAL SPECIFICATIONS

Dimensions

Length	10.7 cm (4.21 in)
--------	-------------------

Width	6.6 cm (2.60 in)
-------	------------------

Height	2.0 cm (0.79 in)
--------	------------------

Case Material

Polybutylene Terephthalate (PBT)

Sample Dish Capacity

7.5 mL, recommended (15 mL, full)

Weight

115 g (4 oz)

Display

64 × 128 graphical

ELECTRICAL AND TIMING CHARACTERISTICS

Operating Temperature

Minimum 4 °C

Typical NA

Maximum 50 °C

Operating Environment

0–90% relative humidity (noncondensing)

Power

2–3 V 16 mm coin cell batteries (life of approximately 3 years)

COMPLIANCE

Manufactured under ISO 9001:2015

EM ISO/IEC 17050:2010 (CE Mark)

3.2 PAWKIT COMPONENTS

Figure 10 shows the features on the top and Figure 11 shows hardware on the bottom of the PAWKIT.

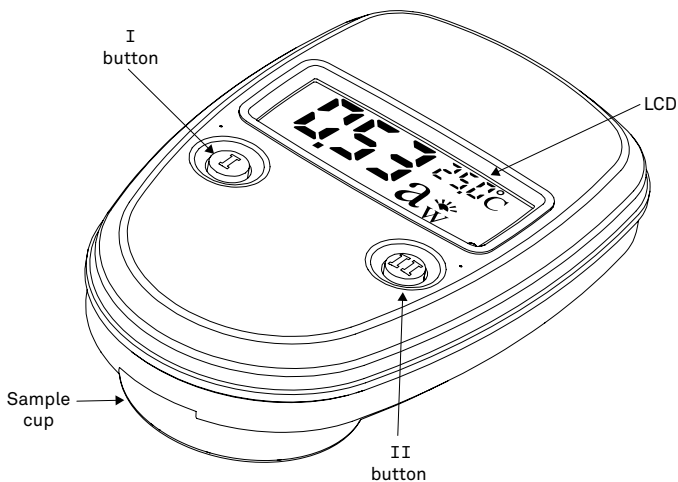


Figure 10 PAWKIT features

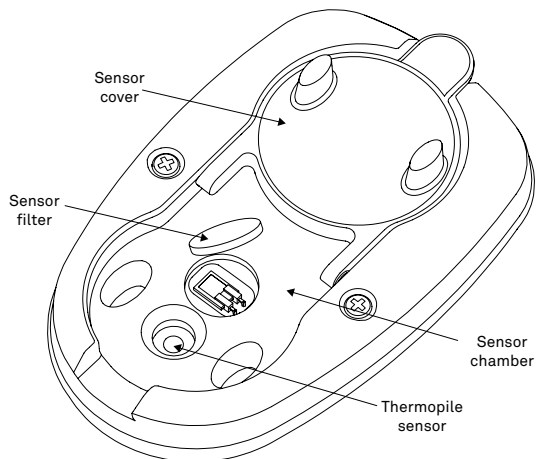


Figure 11 PAWKIT hardware

The PAWKIT uses a capacitance humidity sensor to measure the water activity of a sample. The sensor is suspended behind the sensor filter and uses a special polymer material sandwiched between two microporous electrodes to sense humidity changes. The sensor measures a specific capacitance, which is translated by the software and displayed as water activity on the instrument screen. At equilibrium, the relative humidity of the air in the chamber is the same as the water activity of the sample.

The PAWKIT is accurate to $\pm 0.02 a_w$. For many applications, this accuracy is more than adequate. If a higher accuracy is required, AQUALAB recommends using the AQUALAB 4TE or TDL Water Activity Meter. Contact [Customer Support](#) for more details.

3.3 THEORY

Water is a major component of foods, pharmaceuticals, and cosmetics. Water influences the texture, appearance, taste and spoilage of these products. Quantifying these attributes require two basic types of water analysis: moisture content and water activity (a_w):

3.3.1 MOISTURE CONTENT

Moisture content implies a quantitative analysis to determine the total amount of water present in a sample. There are two primary methods for determining moisture content: loss on drying and Karl Fischer titration. Secondary methods, such as near infrared (NIR) and nuclear magnetic resonance (NMR), may also be used. Moisture content determination is essential in meeting product nutritional labeling regulations, specifying recipes, and monitoring processes.

However, moisture content alone is not a reliable indicator for predicting microbial responses and chemical reactions in materials. The limitations of moisture content measurement are attributed to differences in the intensity that water associates with other components.

3.3.2 WATER ACTIVITY

Water activity (a_w) is a measurement of the energy status of the water in a system. The value indicates how tightly water is bound, structurally or chemically, within a substance. The lower a sample's water activity, the more tightly bound that water is within the sample. The concept of water activity is of particular importance in determining product quality and safety. It predicts safety and stability with respect to microbial growth, chemical and biochemical reaction rates, and physical properties.

Water activity is a measure of the energy status of the water in a system and is a far better indicator of product stability than moisture content. Figure 12 shows how the relative activity of microorganisms, lipids, and enzymes relate to water activity. While other factors, such as nutrient availability and temperature, can affect the relationships, water activity is the best single measure of how water affects these processes.

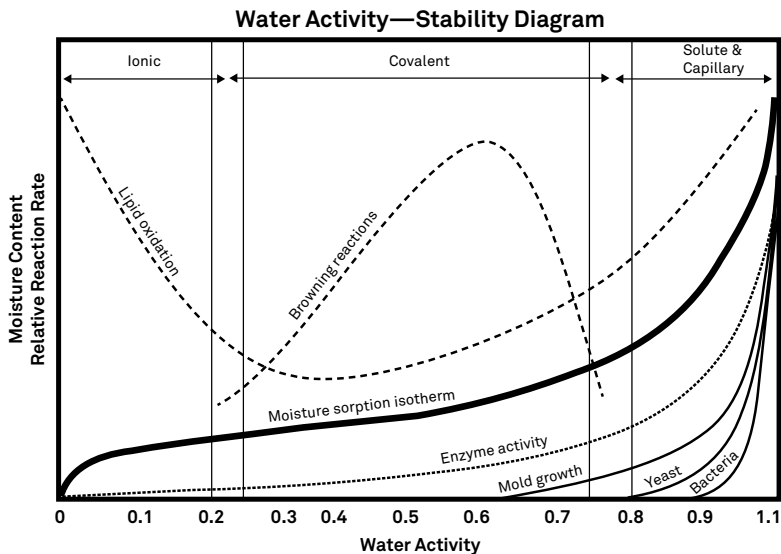


Figure 12 Water activity diagram adapted from Labuza (1976)

Researchers measure the water activity of a system by equilibrating the liquid phase water in the sample with the vapor phase water in the headspace and measuring the relative humidity of the headspace. In the PAWKIT, a sample in a sample cup is placed on a level surface and the PAWKIT is seated on top of the cup, creating a seal. This chamber contains an infrared thermometer and a capacitive humidity sensor. Changes in the electrical capacitance of the polyimide layer of the sensor occur as the relative humidity of the chamber changes. The PAWKIT computes the relative humidity of the headspace by monitoring the change in electrical capacitance. When the water activity of the sample and the relative humidity of the air are in equilibrium, the measurement of the headspace humidity gives the water activity of the sample.

In addition to equilibrium between the liquid phase water and the vapor phase, the internal equilibrium of the sample is important. If a system is not at internal equilibrium, one might measure a steady vapor pressure (over the period of measurement) which is not the true water activity of the system. An example of this might be a baked good or a multicomponent food. Initially out of the oven, a baked good is not at internal equilibrium; the outer surface is at a lower water activity than the center of the baked good. One must wait a period of time in order for the water to migrate and the system to come to internal equilibrium. It is important to keep in mind that by this definition, water activity is an equilibrium measurement.

The water activity of a sample is influenced by factors that affect the binding of water. These factors include temperature, osmotic, matrix, and pressure effects. Typically, water activity is measured at atmospheric pressure, so only the osmotic, temperature, and matrix effects are important.

TEMPERATURE EFFECTS

Temperature plays a critical role in water activity determination. Most critical is the measurement of the difference between sample and the PAWKIT temperature. For water activity measurements to be accurate to 0.02, temperature difference measurements need to be accurate to 0.02 °C. The PAWKIT infrared thermometer measures the temperature of the sample surface. The thermometer is carefully calibrated to minimize temperature errors; however, achieving 0.02 °C accuracy is difficult when temperature differences are large. Most accurate measurements are obtained when the sample is near instrument temperature.

Another effect of temperature on water activity occurs when samples are near saturation. A sample that is close to 1.0 a_w and is warmer than the PAWKIT will cause water to condense within the chamber. This causes errors in the measurement and in subsequent measurements until the condensation disappears. For example, a sample at 0.75 a_w only needs to be 4 °C above the sample chamber temperature to cause condensation to form. For high water activity samples, the operator needs to be aware that condensation can occur if a sample is warmer than the PAWKIT.

The AQUALAB 4TE and TDL have several advantages in having a temperature-controlled chamber. A few major reasons are detailed here.

1. Research purposes.

Researchers can use temperature control to study the effects of temperature on the water activity of a sample, make a comparison of the water activity of different samples independent of temperature, and conduct accelerated shelf-life studies or other water activity studies where temperature control is critical. There are many shelf-life, packaging, and isotherm studies in which temperature control would be very beneficial.

2. Compliance with government or internal regulations for specific products.

Though the water activity of most products varies by less than ± 0.02 per degree Celsius, some regulations require measurement at a specific temperature. The most common specification is 25 °C, though 20 °C is sometimes indicated.

3. Minimization of extreme ambient temperature fluctuations.

If the environmental and PAWKIT temperatures fluctuate by as much as $\pm 5^\circ\text{C}$ daily, water activity readings may vary by $\pm 0.01 a_w$. Temperature control eliminates variations caused by changes in ambient conditions.

The AQUALAB 4TE and TDL models have thermoelectric components installed to allow the instrument to maintain a set chamber temperature.

The water activity of the water in a system is influenced by factors that affect the binding of water. They include osmotic, matrix, and pressure effects. Typically water activity is measured at atmospheric pressure, so only the osmotic and matrix effects are important.

OSMOTIC EFFECTS

Osmotic effects are well known from biology and physical chemistry. Water is diluted when a solute is added. If this diluted water is separated from pure water by a semipermeable membrane, water tends to move from the pure water side through the membrane to the side with the added solute. If sufficient pressure is applied to the solute-water mixture to just stop the flow, this pressure is a measure of the osmotic potential of the solution. Addition of 1 mol of an ideal solute to 1 kg of water produces an osmotic pressure of 22.4 atm. This lowers the water activity of the solution from 1.0 to $0.98 a_w$. For a given amount of solute, increasing the water content of the systems dilutes the solute, decreasing the osmotic pressure, and increasing the water activity. Since microbial cells are high concentrations of solute surrounded by semipermeable membranes, the osmotic effect on the free energy of the water is important for determining microbial water relations and therefore microbe activity.

MATRIX EFFECTS

The sample matrix affects water activity by physically binding water within its structure through adhesive and cohesive forces that hold water in pores and capillaries and to particle surfaces. If cellulose or protein were added to water, the energy status of the water would be reduced. Work would need to be done to extract the water from this matrix. This reduction in energy status of the water is not osmotic, because the cellulose or protein concentrations are far too low to produce any significant dilution of water. The reduction in energy is the result of direct physical binding of water to the cellulose or protein matrix by hydrogen bonding and van der Waal forces. At higher water activity levels, capillary forces and surface tension can also play a role.

3.3.3 WATER POTENTIAL

Water activity is closely related to a thermodynamic property called the water potential, or chemical potential (μ) of water, which is the change in Gibbs free energy (ΔG) when water concentration changes. Equilibrium occurs in a system when μ is the same everywhere in the system. Equilibrium between the liquid and the vapor phases implies that μ is the same in both phases. This means the measured water potential of the vapor phase and can be used to determine the water potential of the liquid phase. Gradients in μ are driving forces

for moisture movement. In an isothermal system, water tends to move from regions of high water potential (high a_w) to regions of low water potential (low a_w). Moisture content is not a driving force for water movement, except in homogeneous materials.

3.3.4 SORPTION ISOTHERMS

Changes in water content affect both the osmotic and matrix binding of water in a product. Therefore a relationship exists between the water activity and moisture content. This relationship is called the sorption isotherm and is unique for each product. Besides being unique to each product, the isotherm changes depending on whether it was obtained by drying or wetting the sample. Typically, large safety margins are built into water content specifications to allow for these uncertainties.

While the sorption isotherm is often used to infer water activity from moisture content, one could easily go the other direction and infer the moisture content from water activity. This is particularly attractive because water activity can be measured much more quickly than water content. This method gives particularly good precision in the center of the isotherm. In order to infer moisture content from water activity, an isotherm for the particular product is needed. AQUALAB can run the isotherm for a fee. Visit aqualab.com/lab-services for more information.

4. SERVICE

This section describes the calibration and maintenance of the PAWKIT. Troubleshooting solutions and customer service information are also provided.

4.1 VERIFICATION AND CALIBRATION

The PAWKIT takes water activity measurements by measuring the change in electrical properties of a special polymer held between two electrodes. Because of the nature of the capacitance humidity sensor, it is important to verify the PAWKIT water activity calibration against known standards to guarantee optimal performance and accuracy. AQUALAB recommends verification daily, once per shift, or before each use. AQUALAB also recommends annual factory calibration to maintain optimal performance.

4.1.1 VERIFICATION STANDARDS

The PAWKIT uses three calibration standards:

- 6.00 mol/kg NaCl ($0.760 a_w$)
- 13.41 mol/kg LiCl ($0.250 a_w$)
- 2.33 mol/kg NaCl ($0.920 a_w$)

Verification standards are specially prepared, unsaturated salt solutions having a specific molality and water activity value that are accurately measurable. The verification standards have been produced under a strict quality assurance regime, and their accuracy is verified by an independent third party instrument. They are very accurate, easy to use, and readily available from AQUALAB. Using verification standards to verify accuracy can greatly reduce preparation errors. For these reasons, AQUALAB recommends using these standards for the most accurate PAWKIT calibration. The verification standards are shelf-stable for 1 year.

To use a verification standard, remove the twist top and pour the contents into an AQUALAB sample cup. Information about the standard value and molality can be found printed on the outside of the plastic vial. If for some reason AQUALAB verification standards are not available, contact [Customer Support](#) for instructions on how to create a verification standard salt solution.

The capacitance sensor can exhibit hysteresis. When a high water activity standard or sample is measured, this may influence future readings. After measuring a high water activity sample, allow 1 h for the sensor to dry. Additionally, desiccant or activated carbon can be used to accelerate the drying process.

NOTE: To avoid inaccurate water activity readings, verification standards should be used once immediately after opening and not stored in sample cups for repeated use.

4.1.2 VERIFY CALIBRATION

Use the following steps to verify the PAWKIT calibration.

NOTE: An error code of 9.99 at any time during the process indicates that the sensor has failed and that the instrument needs to be serviced. Refer to [Section 4.4](#) for shipping instructions.

1. Using a vial of the 0.760 a_w NaCl standard, empty the entire contents of the vial into a sample cup.
2. Place the PAWKIT over the sample cup as described in the [Section 2.1.2](#).
3. Press the I button to take a reading.

If it is reading the correct water activity (± 0.02), the PAWKIT needs no adjustment for this standard. Skip to [step 7](#).

If the first reading was not the correct water activity (± 0.02), clean the PAWKIT according to the instructions in [Section 4.2](#) and take a second reading.

If the second reading is the correct water activity (± 0.02), the PAWKIT needs no adjustment at this time. Skip to [step 7](#). If the second reading is not correct, continue to the next step.

4. Press the II button (on the right) once to get the **Calibration Mode** screen ([Figure 13](#)).

Once the reading is finished, the II button (on the right) activates. The II button is only active until the PAWKIT shuts itself off.

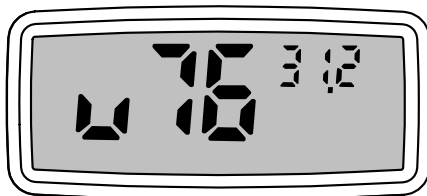


Figure 13 Calibration Mode screen

[Figure 13](#) shows that the PAWKIT is ready to have its calibration adjusted up for the 0.76 standard. The number in the upper right corner indicates the water activity measurement that the PAWKIT just read. Press the II button to scroll through the other selections. They are: u76, d76, u25, d25, Sto, u92, and d92. The u and d before each number stand for **up** or **down** adjustment for each standard. The numbers (25, 76, and 92) correspond to the water activity of a verification standard (0.76, 0.25, and 0.92 a_w). The Sto position stores a reading.

EXAMPLE:

If the NaCl reading is lower than it should be, press the II button to scroll to u76 (adjust up for 0.76 standard). If the reading is higher than it should be, scroll to d76 (adjust down for 0.76 standard).

NOTE: Scroll feature is continuous, so if desired adjustment screen is passed, keep pressing the II button until cycled back to the correct screen.

5. Adjust the calibration.
 - a. Scroll to the proper screen for calibration adjustment.
 - b. Press the **I** button to adjust the correct value.
Each time the **I** button is pressed, the value in the corner changes by an increment of 0.01.
6. Store the new calibration value.
 - a. Scroll to the proper screen for calibration adjustment.
 - b. Press the **II** button to scroll until **sto** appears in the lower right corner.
 - c. Press the **I** button to store the new calibration value.
 - d. Return to the main screen and begin a new measurement.

NOTE: If **sto** is not pressed, no change is made to the PAWKIT calibration.

7. Verify with a second standard, either the 0.25 or the 0.92 standard.
Choose the standard that is closest to the water activity range of the sample material being tested. In other words, if it is normally higher than $0.76 a_w$, use the 0.92 standard. If it is normally lower than $0.76 a_w$, use the 0.25 standard. If the PAWKIT measures the second standard correctly (± 0.02), begin testing the product. If it does not measure correctly, repeat [step 3](#) through [step 7](#) for the second standard.

8. If the calibration routine is inadvertently entered, keep pressing the **II** button until the main screen appears.

NOTE: The 0.76 standard adjustment adjusts the calibration intercept, while the 0.25 and 0.92 adjusts the slope. Changes in the intercept are more likely to occur than changes in the slope, so the 0.76 verification check is the most important and should be done more frequently.

Review the graphical representation of the calibration routine in [Figure 14](#).

SERVICE

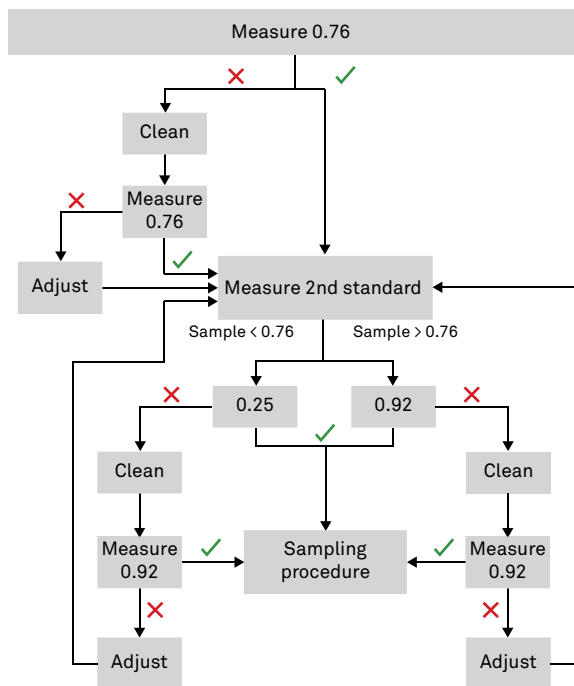


Figure 14 Calibration routine flowchart

4.2 MAINTENANCE

The PAWKIT may be returned to AQUALAB for maintenance in the following areas: system inspection, parts replacement, and instrument cleaning. Replacement parts can also be ordered from AQUALAB. Contact Customer Support for more information.

4.2.1 CLEANING

The accuracy of the PAWKIT relies on keeping the instrument clean. Dust and sample debris can contaminate the sample chamber and must be removed regularly. Follow the instructions in this section carefully to clean the instrument.

Materials Needed for Cleaning

Before beginning the cleaning process, prepare the following items:

- 3 or more cleaning swabs
- Deionized water
- IPA cleaning solution

IMPORTANT:

- DO NOT get the capacitance humidity sensor filter wet or touch it with your fingers as water and oils will slow the sensor response and increase the reading duration and may cause the sensor to read incorrectly. If sample material spills on the sensor filter, it must be replaced following the instructions in Section 4.2.2. Preventing contamination of this filter is critical, as it will influence the measurement
- Wash your hands or wear gloves before starting the cleaning process.

Cleaning the Outer Enclosure

- **LCD Screen:** Use only a soft cotton cloth to clean the LCD. Do not use tissues, as they may scratch the surface.
- **Outer Case:** Clean the rest of the outer enclosure with a soft cotton cloth.

Cleaning the Sensor Chamber

1. Wash

- Apply two drops of deionized water to a fresh cleaning swab.
- Use the swab to clean the thermopile sensor lens, the chamber area, and recess of the sample chamber (where the rim of the cup goes).
- Use additional swabs if necessary to ensure thorough cleaning.

2. Clean

- Apply two drops of IPA cleaning solution to a fresh swab.
- Clean the thermopile sensor lens, chamber area, and recess of the sample chamber (where the rim of the cup goes).
- Ensure all surfaces are free of dirt, lint, and particles that could cause contamination. Repeat with fresh swabs if necessary.

3. Dry

- Use dry, fresh swabs to gently and thoroughly dry the PAWKIT.

4. Visual Inspection

- Inspect the instrument for cleanliness. If the chamber remains dirty or if the drying swab shows discoloration or debris, repeat the cleaning steps.

By following these steps, you will maintain the PAWKIT's accuracy and ensure consistent performance.

NOTE: Do NOT reuse swabs.

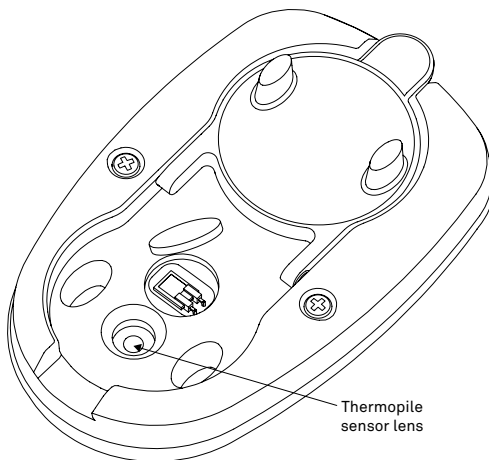


Figure 15 Thermopile sensor lens

4.2.2 FILTER REPLACEMENT

The porous white humidity sensor filter may periodically need to be replaced if it becomes dirty. To purchase replacement sensor filters, please visit store.aqualab.com

1. To remove the sensor filter use a knife or needle point to gently pry up the edge of the filter (Figure 16).
2. Be careful of the capacitance humidity sensor (Figure 17) when placing the new filter. Use gloves to handle the filters as finger oils can be transferred to the filter and contaminate the filter.
3. Place one edge of the filter into the sensor chamber without touching the sensor. Use a finger to push the filter into place.

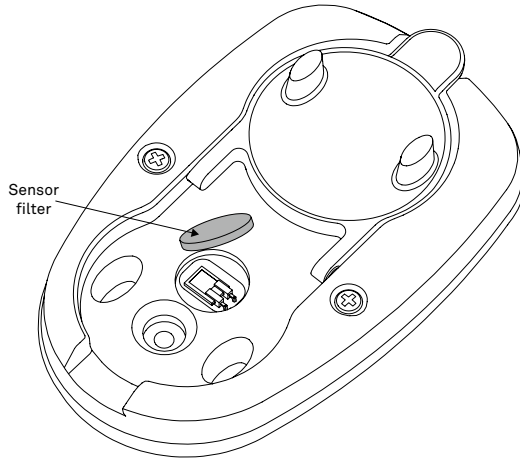


Figure 16 Filter removal

⚠ CAUTION

DO NOT touch the capacitance humidity sensor because it is extremely fragile.

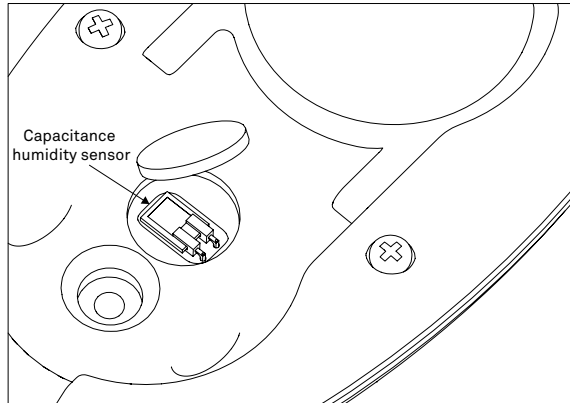


Figure 17 Capacitance humidity sensor

4.2.3 BATTERY REPLACEMENT

The PAWKIT uses two lithium-ion battery cells that should last for several years. If the battery charge is low, a low-battery indicator icon (Figure 18) will persistently appear in the lower right corner of the screen.

NOTE: If the low battery indication appears periodically and is not persistent, it does not necessarily mean the battery needs to be replaced.

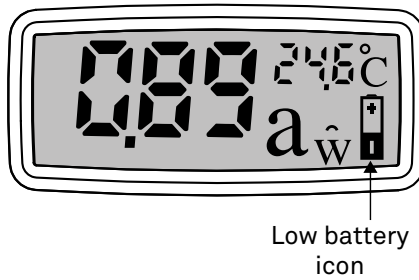


Figure 18 Low battery icon

Follow steps 1 through 6 below to replace battery.

1. Remove the PAWKIT bottom by unscrewing the two screws.

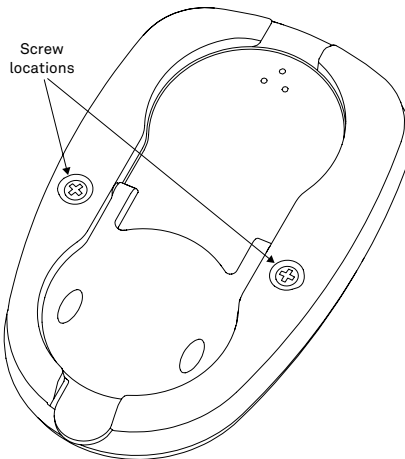


Figure 19 Battery compartment screws

2. Separate the stainless steel top and elastomer (which contains the batteries) from the white plastic bottom.

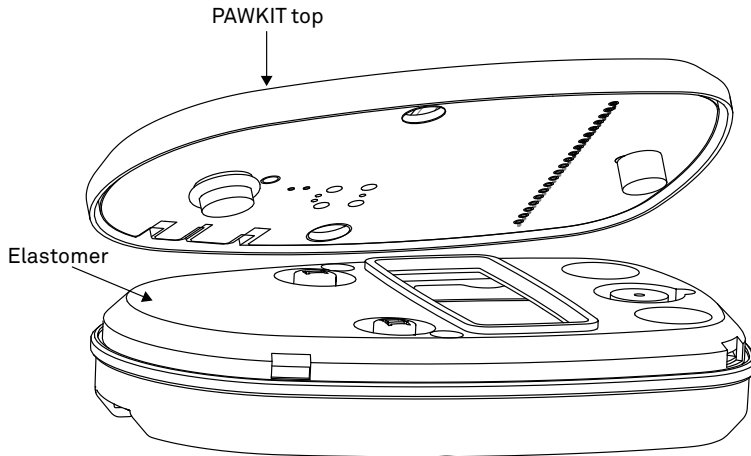


Figure 20 Separate top and elastomer

3. Remove the old lithium-ion batteries.
4. Replace with new CR1632 or equivalent 3 V lithium coin cells.

Make sure to orient the batteries so positive (+) contact is facing down into the elastomer pocket. Make sure the two small springs (Figure 21) which make contact between the (+) battery terminal and the circuit board are in place.

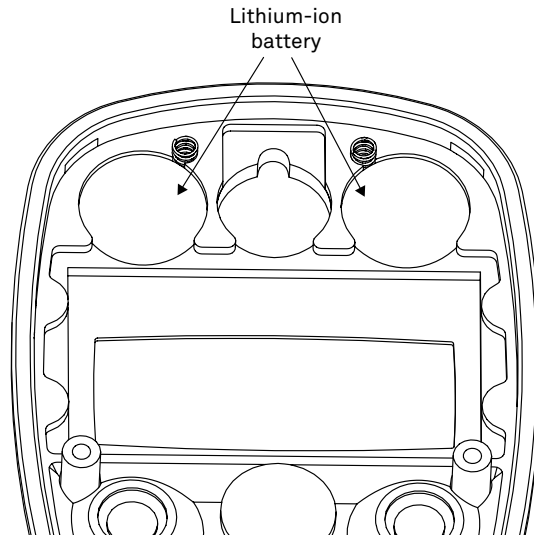


Figure 21 Lithium-ion batteries and contact springs

5. Replace the circuit board/plastic bottom in the elastomer top.
6. Tighten the two screws to complete the assembly of the PAWKIT.

4.3 REPAIRS

AQUALAB repairs manufacturer defects and instruments within the 1-year warranty at no charge. Repairs outside of the warranty window are charged based on cost of parts, labor, and shipping. An extra fee may be charged for rush work. Contact [Customer Support](#) for an estimated repair cost. AQUALAB has loaner instruments available for a fee while the PAWKIT is being serviced.

All PAWKIT units returning to AQUALAB for servicing must be accompanied with a Return Merchandise Authorization (RMA) number. Prior to shipping the instrument, contact Customer Support to obtain an RMA number. The following steps should help to ensure the safe shipping and processing of the PAWKIT:

1. Ship the PAWKIT in its original hard carry case with suspension packaging.
If it is not possible, use a box with at least 4 in of space between the instrument and each wall of the box.
2. Place the PAWKIT in a plastic bag to avoid disfiguring marks from the packaging.
3. If the original packaging is not available, pack the box moderately tight with packing material (e.g., Styrofoam® peanuts or bubble wrap), ensuring the instrument is suspended in the packing material.
4. On the RMA form, please verify the ship to and bill to information, contact name, and problem description. If anything is incorrect, please contact [Customer Support](#).
5. Tape the box in both directions for added support.
6. Include the RMA number in the attention line on the shipping label.

A Certificate of Calibration will be issued upon completion of the work.

4.4 TROUBLESHOOTING

[Table 2](#) lists common problems and their solutions. If the problem is not listed or these solutions do not solve the issue, contact [Customer Support](#).

Table 2 Troubleshooting the PAWKIT

Problem	Possible Solutions
Error Code 9.99	The sensor has failed. Refer to Section 4.4 for further instructions on how to return the PAWKIT for repair.
Constantly offsetting or consistently high readings.	<p>The instrument may be dirty. Refer to Section 4.2.1 for directions on cleaning.</p> <p>The filter may be dirty. Refer to Section 4.2.2 for directions on changing the filter.</p>

Table 3 Troubleshooting the PAWKIT (continued)

Problem	Possible Solutions
Visible debris or staining	Refer to Section 4.2.1 for directions on cleaning and Section 4.2.2 for directions on changing the filter. Refer to Section 2.1.1 on sample preparation. Be careful to not get sample on the rim of the cup. Wipe the rim before placing the instrument on the cup.
Corrosion on thermopile	Contact Customer Support for repair. Refer to Section 4.4 for detailed instructions.
Instrument is unstable when depressing a button	Move the instrument to a flat surface.

4.5 CUSTOMER SUPPORT

NORTH AMERICA

Customer service representatives are available for questions, problems, or feedback Monday through Friday, 7 am–5 pm Pacific time.

Email: support@aqualab.com
sales@aqualab.com

Phone: +1.509.332.5601

Fax: +1.509.332.5158

Website: aqualab.com

EUROPE

Customer service representatives are available for questions, problems, or feedback Monday through Friday, 8:00 to 17:00 Central European time.

Email: support.europe@aqualab.com
sales.europe@aqualab.com

Phone: +49 89 208049202

Website: aqualab-eu.de

If contacting AQUALAB by email, please include the following information:

SERVICE

Name	Email address
Address	Instrument serial number
Phone	Description of the problem

NOTE: For products purchased through a distributor, please contact the distributor directly for assistance.

4.6 TERMS AND CONDITIONS

By using AQUALAB instruments and documentation, you agree to abide by the AQUALAB by Addium, Inc. USA Terms and Conditions. Please refer to addium.io/en/company/terms-and-conditions for details.

REFERENCE

Labuza, Ted P., K. Acott, S. R. TatiNI, R. Y. Lee, Jv Flink, and W. McCall. "Water activity determination: a collaborative study of different methods." *Journal of Food Science* 41, no. 4 (1976): 910-917.

INDEX

A

accuracy 9
AQUALAB 11

B

battery
 lithium-ion 22
 replacement 22
beeper 2, 4
binding 13–15
buttons
 I button 6
 adjust calibration 18
 restart measurement 6
 verification, start reading 17
 wakes up PAWKIT 8
 II button 17–18
 calibration mode screen 17
 scroll to desired calibration mode 17
 turn on PAWKIT 6

C

calibration 16–18
 precautions 2
 standards (LiCl, NaCl) 16–18
 d25 17
 d76 17
 d92 17
 molality 16
 Sto 17
 u25 17
 u76 17
 u92 17
capillaries 14
cleaning. *See* maintenance, cleaning
close sample chamber lid. *See* sample
 chamber lid, close
cold sample. *See* temperature

compliance 10
 CE 10
cosmetics. *See* theory
customer support 26

D

d25. *See* calibration, standards
d76. *See* calibration, standards
d92. *See* calibration, standards
display
 cleaning LCD 20–21
 LCD 6

E

electrical and timing characteristics 10
 operating environment 10
 operating temperature 10
 power requirement 10
enzymes. *See* theory; water activity
equilibrium. *See* theory; water activity
 water activity and relative humidity 13

F

features. *See* PAWKIT, features
filter 21–22
 cleaning 20–21
 replacement 21–22

H

hardware. *See* PAWKIT, hardware
humidity
 related to water activity 11

L

LiCl. *See* calibration, standards
lipids. *See* water activity, lipids

liquid phase water.

See water activity, liquid phase water

lithium-ion batter.

See maintenance, battery replacement

M

maintenance

battery replacement 22

cleaning 19

lithium-ion battery 22

repairs. See maintenance, repairs

measurement

taking 4

time 7

moisture content 11

definition 11

loss on drying 11

quantitative analysis 11

versus water activity 12

molality of standards.

See calibration, standards, molality

N

NaCl. See calibration, standards

O

OFF

turn PAWKIT OFF 8

ON

turn PAWKIT ON 6

open sample chamber lid.

See sample chamber lid, open

P

PAWKIT

accessories 1

features 10

hardware 11

OFF 8

ON 6

operation 2–8

precautions

capacitance humidity sensor 22–23

sampling 2

preparation

for operation 2

samples 3

pressure effect.

See water activity, factors in determining

R

references

relative humidity

affects cleaning 20

features 11

operating environment 10

repairs. See maintenance, repairs

S

sample chamber lid

close 8

open 4

sample cups

filling level 3

lids 3

stainless steel 11

sampling

environment 3

inserting sample 5–6

location 2

multicomponent 3, 13

preparation 3

saturated sample.

See temperature, wet sample, cautions

sensor, capacitance humidity 11

damage 22

filter

cleaning 21

verification 16

sorption isotherm 15

sorption isotherms. *See also* theory
 specifications 9–10
 measurement specifications 9
 accuracy 9
 read time 9
 sample temperature 9
 water activity 9
 physical specifications 9
 case material 9
 dimensions 9
 sample dish capacity 9
 weight 9
 spoilage. *See* theory, perishability
 stability diagram.
 See theory, stability diagram

T

temperature
 cold sample 4
 effects on water activity 12
 equilibrium 11
 hot sample 3–4
 wet sample, cautions 15
 terms and conditions 27
 theory
 binding 13–15
 capillaries 14
 cosmetics 11
 enzymes 12
 equilibrium
 water activity and relative humidity 11
 factors in determining
 pressure effects 13
 Gibbs free energy 15
 homogenous 3
 humidity 11–12
 lipids 12
 liquid phase water 12
 loss on drying 11
 microbial growth 12
 multicomponent food 13
 osmotic effects 14
 perishability 11

 pharmaceuticals 11
 quantitative analysis 11
 regulations 11–13
 thermodynamic 15
 water activity 12–14
 thermodynamic. *See* water potential
 time. *See* measurement, time
 troubleshooting
 corrosion on thermopile 25
 error code 9.99 25
 high readings 25
 offsetting 25
 unstable when depressing button 25
 visible debris or staining 25

W

water activity 12–14
 effect on food 11–13
 enzymes 12
 equilibrium 11
 lipids 12
 liquid phase water 12
 microbial growth 12
 multicomponent food 13
 osmotic effects 14
 perishability 11. *See also* theory
 regulations 11–13
 stability diagram 12
 water potential
 Gibbs free energy 15
 homogeneous 3
 thermodynamic property 15